

632.95

261

H

ISO

EPA-600/8-80-038
June 1980

ANALYSIS OF PESTICIDE RESIDUES IN HUMAN AND ENVIRONMENTAL SAMPLES

**A COMPILATION OF METHODS
SELECTED FOR USE IN
PESTICIDE MONITORING PROGRAMS**

REVISIONS BY

THE ASSOCIATION OF OFFICIAL ANALYTICAL CHEMISTS

JOSEPH SHERMA
MORTON BEROZA



111962

CONTRACT NO. 68-02-2474

EDITED BY

RANDALL R. WATTS
REVISED DECEMBER 1979

EDITORIAL PANEL

R.G. LEWIS
R.F. MOSEMAN
D.W. HODGSON

**U.S. ENVIRONMENTAL PROTECTION AGENCY
HEALTH EFFECTS RESEARCH LABORATORY
ENVIRONMENTAL TOXICOLOGY DIVISION
RESEARCH TRIANGLE PARK, NORTH CAROLINA**

TABLE OF CONTENTS*

<u>Section</u>	<u>Subject</u>
1.	Introduction
2.	Collection, preservation, and storage of samples
	I. General comments
	II. Sample containers
	III. Storage of samples
	IV. Sampling, general
3.	Miscellaneous information
	A. Cleaning of laboratory glassware
	B. Preparation, storage, and use of pesticide analytical standards
	C. General purity tests for solvents and reagents
	D. Evaluation of quality of Florisil
	E. Limits of detectability - blood and adipose tissue
4.	Gas-liquid chromatography
	A. Electron capture detection
	(1) Description of instrument and accessories
	(2) Columns
	(3) Detector
	(4) Chromatography of sample
	(5) Quantitation and interpretation
	(6) Tables and figures
	(7) Support-bonded Carbowax 20M columns

*Analytical methods designated by a star have been subjected to interlaboratory study.

4. (Cont.)

B. Flame photometric detection

- (1) Description of component modules
- (2) Columns
- (3) Detector
- (4) Sample quantitation and interpretation
- (5) Tables and figures

C. Hall electrolytic conductivity detector

- (1) Description of instrument
- (2) Columns
- (3) Detector
- (4) Sample quantitation and interpretation
- (5) Retention data and chromatograms of carbamates on Carbowax-modified supports

D. Nitrogen-phosphorus (N-P) detector

5. Chlorinated hydrocarbon pesticides and metabolites

A. In human tissues and excreta

(1) Adipose tissue

★ (a) Modification of Mills, Onley, Gaither method for multiple chlorinated pesticides and metabolites

★ (b) Hexachlorobenzene (HCB) and mirex, and HCB confirmation

(2) Micro methods

(a) Liver, kidney, bone, marrow, adrenal, gonads

(b) Brain and human milk

- 5. (Cont.)
 - (3) Blood
 - ★ (a) Multiple chlorinated pesticides
 - (b) Pentachlorophenol
 - (4) Urine
 - (a) Pentachlorophenol and other chlorinated phenols
 - (b) DDA in urine
 - (c) 2,4-D and 2,4,5-T
 - (5) Human blood and environmental samples
 - (a) Kepone
 - B. Cleanup by gel permeation chromatography (GPC)
- 6. Organophosphorus pesticides and metabolites
 - A. In human tissues and excreta
 - (1) General comments
 - (2) Urine
 - (a) Determination of metabolites or hydrolysis products of organophosphorus pesticides
 - (b) Para-nitrophenol
 - (3) Blood
 - ★ (a) Cholinesterase activity
- 7. Carbamate pesticides and metabolites
 - A. 1-Naphthol in urine
- 8. Sampling and analysis of air for pesticides
 - A. Sampling
 - B. Analysis

9. Analysis of nonpesticide pollutants: PCBs and TCDD
 - A. Introduction to the analysis of PCBs
 - B. Determination of PCBs in human milk
 - (1) Macro method
 - (2) Micro methods
 - C. Separation of PCBs from organochlorine pesticides
 - D. TLC semi-quantitative estimation of PCBs in adipose tissue
 - E. Typical PCB chromatograms
 - F. Tables of relative retention and response ratios of PCBs
 - G. Analysis of TCDD in milk, liver, fish, water, and sediment
10. Analysis of water
 - A. Sampling and analysis of water for pesticides
 - B. Determination of some free herbicides in water
11. The analysis of soils, housedust, and sediment
 - A. Organochlorine insecticides in soils and housedust
 - B. Organochlorine and organophosphorus insecticides in bottom sediment
 - C. Carbamate pesticides in soil
12. Confirmatory procedures
 - A. Confirmation and determination of organochlorine pesticides in human tissue and milk
 - B. Thin-layer chromatography
 - C. p-Values

- D. Derivatization techniques
 - (1) Microscale alkali treatment for confirmation of sample cleanup
 - (2) Perchlorination of PCBs
 - E. Infrared spectroscopy
 - F. Polarography
13. High performance liquid chromatography

APPENDIX

<u>Section</u>	<u>Subject</u>
I.	Maintenance and repair of instruments
II.	A. Analytical quality control
VI.	Block diagram of tentative tissue, excreta and method selection for abnormal pesticide exposure cases
VII.	Pesticide analytical reference standards repository
VIII.	Future manual revisions - very important